

Success

Column Installation Steps

1. Check traps, carrier gas, septum, liner
2. Place the nut and ferrule on the column and carefully cut the column end
3. Install the column into the injector
4. Turn on the carrier gas
5. Install the column into the detector
6. Inspect for leaks
7. Confirm carrier gas flow and proper column installation
8. Condition the column
9. Accurately set the carrier gas velocity
10. Bleed test
11. Run test mix

J&W Technical Support: 800-552-0413
916-985-7888

Web Site: <http://www.jandw.com>

